

INTRODUCTION

NARSIPUR is a renowned ISO 9001:2015 certified company engaged in the manufacture and sale of various hygiene and sanitation chemicals and products. ENCIDAY has been developed by NARSIPUR as a new generation Milk analyser for daily / alkaline cleaner with the objective of easily and effectively removing all types of deposits.

ADVANTAGES

- Fast action with very long residual activity
- Provides superior cleaning action
- Non staining
- Highly stable
- Almost completely bio degradable
- Low corrosivity
- Safe and ecofriendly

USAGE for milk analyser cleaning

Use dilution in Distilled Water for cleaning the automatic Milk analyser.
Daily 2% (1:50) solution

PREPARATION of diluted working solution

Use pipette to add 10 ml ENCIDAY slowly to a beaker or suitable glassware with 490 ml of Distilled Water. Stir and pour into labelled containers.

CAUTION for use:

Keep the concentrate out of reach of children.

If the concentrate is accidentally spilled on the skin, flush the area immediately with running water.

Never transfer the concentrate into unlabelled or improperly labelled containers.

PRESENTATION

In plastic bottles of 200 ml each
